

目录

乙酰唑胺-Acetazolamide.....	3
L-抗坏血酸钠-Sodium Ascorbate.....	3
L-抗坏血酸-Ascorbic acid.....	3
二丙酸倍他米松-Betamethasone dipropionate.....	3
盐酸溴己新-Bromhexine hydrochloride.....	4
盐酸度硫平-Dosulepin hydrochloride.....	4
茶苯海明-Dimenhydrinate.....	4
盐酸多塞平-Doxepin Hydrochloride.....	4
盐酸乙胺丁醇-Ethambutol Hydrochloride.....	5
碘克沙酸-Ioxaglic acid.....	5
醋酸甲羟孕酮-Medroxyprogesterone acetate.....	5
来酸甲麦角新碱-Methylethergometrine Maleate.....	5
诺孕酯-Norgestimate.....	6
福尔可定-Pholcodine.....	6
癸酸睾酮-Testosterone Decanoate.....	6
异己酸睾酮-Testosterone Isocaproate.....	6
庚酸睾酮-Testosterone enantate.....	7
培美曲塞二钠七水合物-Pemetrexed disodium heptahydrate.....	7
利塞膦酸钠 2.5 水合物-Risedronate sodium 2.5-hydrate.....	7
硫酸庆大霉素- Gentamicin sulfate.....	7
硫酸氢氯吡格雷-Clopidogrel hydrogen sulfate.....	8
吗多明-Molsidomine.....	8
度他雄胺-Dutasteride.....	8
米诺地尔-Minoxidil.....	8
2,4-二氯苄醇-2,4-Dichlorobenzyl alcohol.....	9
卡比多巴-Carbidopa.....	9
异麦芽酮糖醇-Isomalt.....	9
二水合盐酸纳洛酮-Naloxone hydrochloride dihydrate.....	9
磺胺二甲氧嗪-Sulfadimethoxine.....	10
氯己定二葡萄糖酸盐溶液- Chlorhexidine digluconate solution.....	10
甲基苯巴比妥-Methylphenobarbital.....	10
磺胺嘧啶-Sulfadiazine.....	10
甲苯磺丁脲-Tolbutamide.....	11
尼泊金甲酯-Methyl parahydroxybenzoate.....	11
尼泊金丙酯-Propyl parahydroxybenzoate.....	11
烟酸-Nicotinic acid.....	11
盐酸布比卡因-Bupivacaine hydrochloride.....	12
尼泊金丁酯-Butyl parahydroxybenzoate.....	12


尼泊金乙酯-Ethyl parahydroxybenzoate.....	12
盐酸氟奋乃静-Fluphenazine dihydrochloride.....	12
尼泊金乙酯钠盐- Sodium ethyl parahydroxybenzoate.....	13
尼泊金甲酯钠盐-Sodium methyl parahydroxybenzoate.....	13
尼泊金丙酯钠盐-Sodium propyl parahydroxybenzoate.....	13
醋酸氟卡胺-Flecainide acetate.....	13
左卡尼汀- Levocarnitine.....	14
比卡鲁胺-Bicalutamide.....	14
氢溴酸西酞普兰-Citalopram hydrobromide.....	14
盐酸多佐胺-Dorzolamide hydrochloride.....	14
奥卡西平-Oxcarbazepine.....	15
盐酸溴己新-Bromhexine hydrochloride.....	15
苯巴比妥钠-Phenobarbital sodium.....	15
钆布醇一水合物-Gadobutrol monohydrate.....	15
恩替卡韦一水合物-Entecavir monohydrate.....	16
草酸艾司西酞普兰-Escitalopram oxalate.....	16
他卡西醇一水合物- Tacalcitol monohydrate.....	16
甲酸-Formic acid.....	16
他克莫司一水合物-Tacrolimus monohydrate.....	17
艾司西酞普兰-Escitalopram.....	17
替莫唑胺-Temozolomide.....	17
蔗糖-Sucrose, liquid.....	17
盐酸尼卡地平-Nicardipine hydrochloride.....	18
替考拉宁-Teicoplanin.....	18
富马酸卢帕他定-Rupatadine fumarate.....	18
替加环素-Tigecycline.....	18
依那西普-Ethosuximide.....	19
唑来磷酸一水化合物-Zoledronic acid monohydrate.....	19
骨化二醇一水合物-Calcifediol monohydrate.....	19
盐酸氟哌噻吨-Flupentixol dihydrochloride.....	19
硫酸阿扎那韦-Atazanavir sulfate.....	20
盐酸多奈哌齐-Donepezil hydrochloride.....	20
盐酸多奈哌齐一水合物-Donepezil hydrochloride monohydrate.....	20
巯嘌呤一水合物-Mercaptopurine monohydrate.....	20
硫酸鱼精蛋白-Protamine sulfate.....	21

乙酰唑胺-Acetazolamide

有关物质: Synergi Polar-RP (4.6x150,4μm)

EUROPEAN PHARMACOPOEIA 10

Acetazolamide

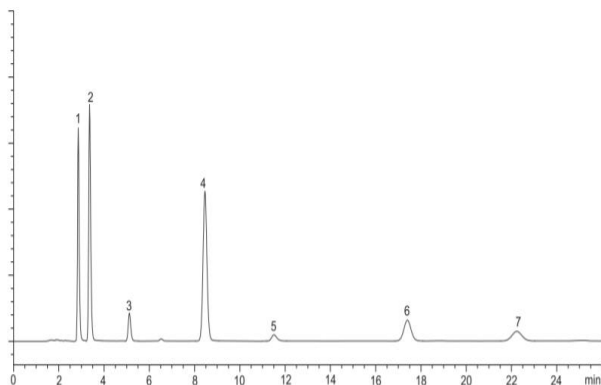


1. impurity E 3. impurity B 5. impurity C 7. impurity F
2. impurity D 4. acetazolamide 6. impurity A

L-抗坏血酸钠-Sodium Ascorbate

有关物质: Sphereclone NH2 (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Sodium ascorbate

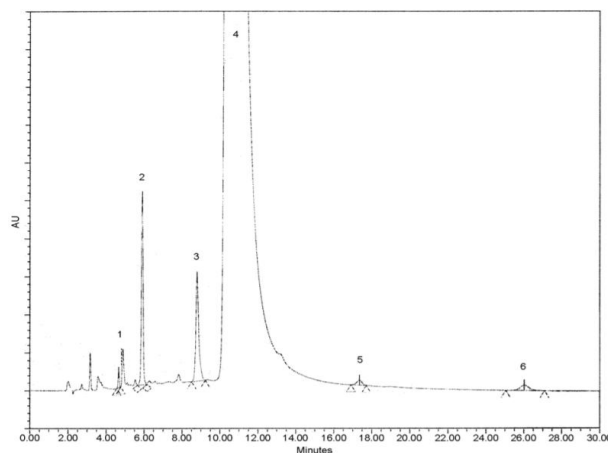


1. impurity D 2. impurity H 3. impurity F 4. ascorbic acid 5. impurity C 6. impurity G

L-抗坏血酸-Ascorbic acid

有关物质: Sphereclone NH2 (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Ascorbic acid

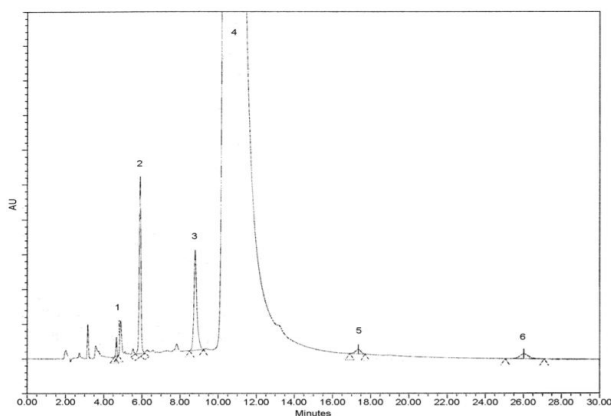


1. impurity D 2. impurity H 3. impurity F 4. ascorbic acid 5. impurity C 6. impurity G

二丙酸倍他米松-Betamethasone dipropionate

有关物质和含量: Luna C18(2)(2.0x100,2.5μm)

EUROPEAN PHARMACOPOEIA 10

Betamethasone dipropionate

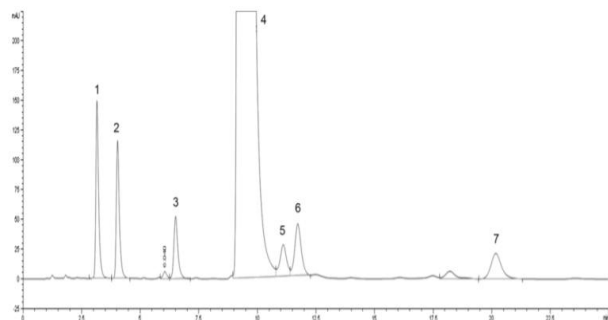


1. impurity B 3. impurity D 5. impurity I 7. impurity G
2. impurity C 4. betamethasone dipropionate 6. impurity E

盐酸溴己新-Bromhexine

hydrochloride


有关物质和含量: Kinetex C18 (2.1x150,2.6μm)

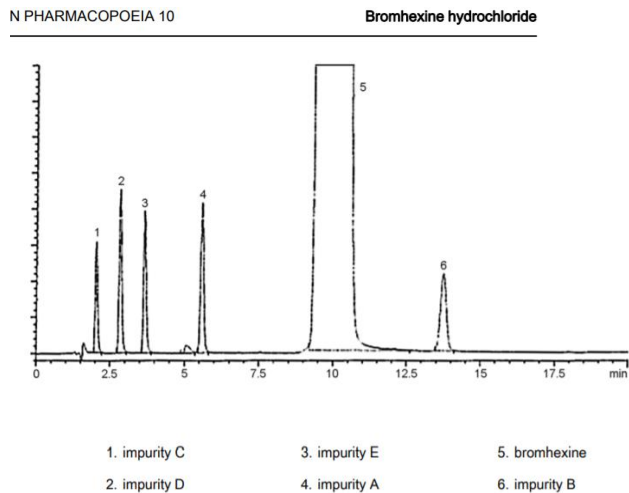

盐酸度硫平-Dosulepin


hydrochloride

有关物质: Luna CN (4.6 x250,5μm)

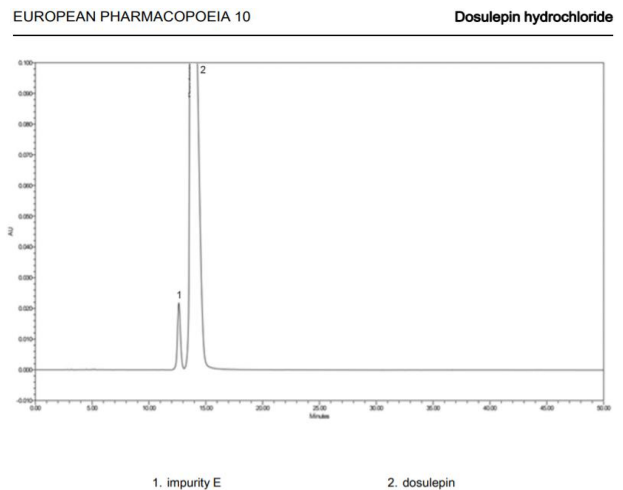


茶苯海明-Dimenhydrinate

有关物质: Luna C18 (4.6x250,5μm)

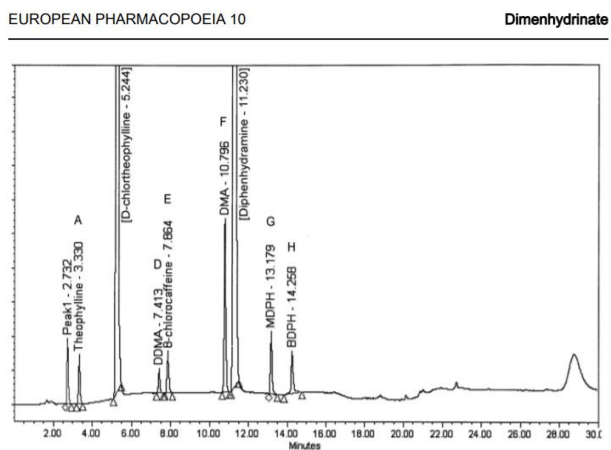

盐酸多塞平-Doxepin Hydrochloride

有关物质: Luna C18 (4.6 x250,5μm)

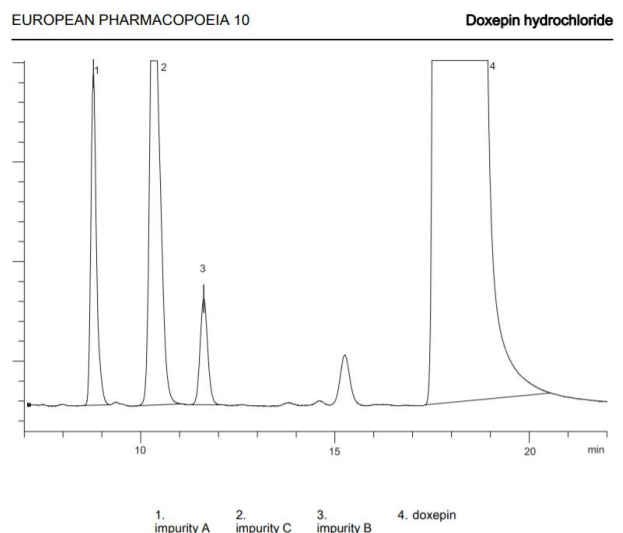


盐酸乙胺丁醇-Ethambutol

Hydrochloride

有关物质: Luna C18(2) (4.6 x100,3μm)

EUROPEAN PHARMACOPOEIA 10

Ethambutol hydrochloride

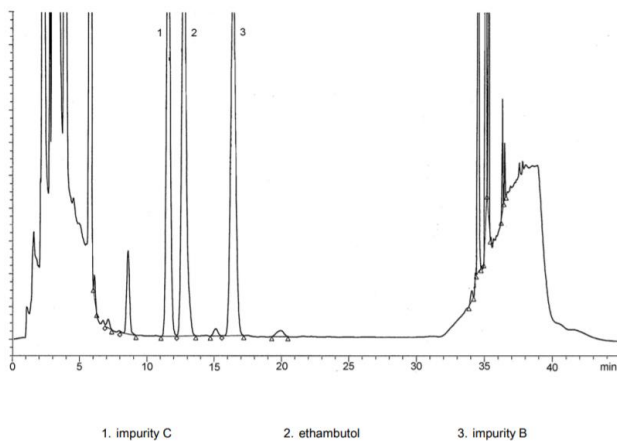


碘克沙酸-Ioxaglic acid

有关物质: Luna C8 (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Ioxaglic acid

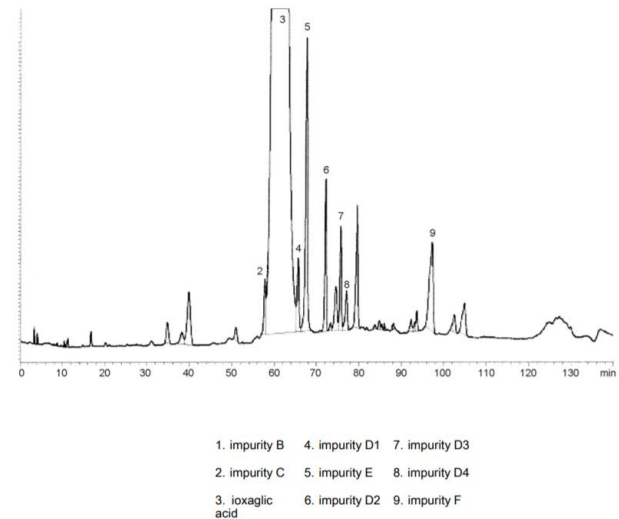

醋酸甲羟孕酮


-Medroxyprogesterone acetate

有关物质: Prodigy ODS-3 (3.0 x250,5μm)

杂质 F: Luna C18 (4.6 x100,3μm)

EUROPEAN PHARMACOPOEIA 10

Medroxyprogesterone acetate

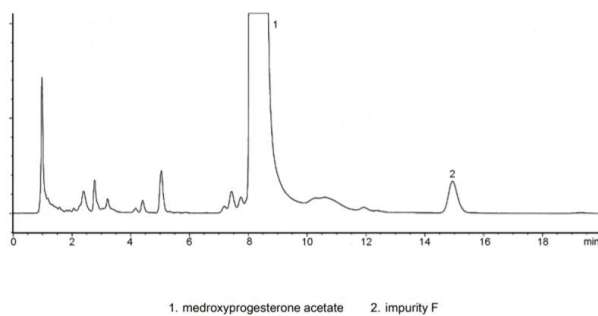


来酸甲麦角新碱

-Methylethergometrine Maleate

有关物质: Ultracarb 5 ODS (30) (4.6 x100,5μm)

EUROPEAN PHARMACOPOEIA 10

Methylethergometrine maleate

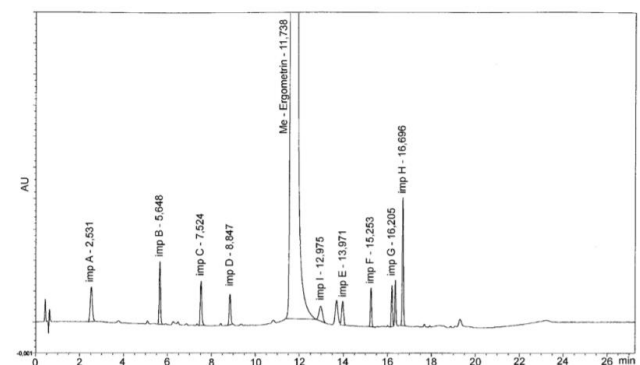


诺孕酯-Norgestimate

有关物质: Luna C18 (4.6 x100,5µm)

EUROPEAN PHARMACOPOEIA 10

Norgestimate

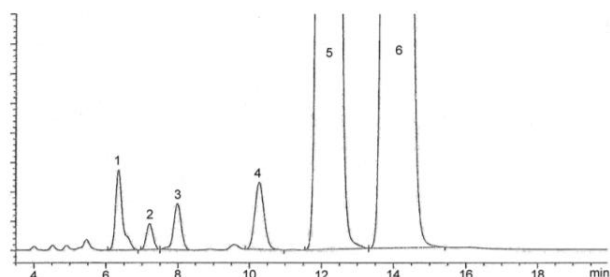


- | | | |
|---------------|---------------|-------------------------------|
| 1. impurity B | 3. impurity D | 5. (Z)-isomer of norgestimate |
| 2. impurity C | 4. impurity A | 6. (E)-isomer of norgestimate |

福尔可定-Pholcodine

有关物质: Luna Phenyl-Hexyl (4.6 x75,3µm)
或 Kinetex Phenyl-Hexyl (4.6 x75,2.6µm)

EUROPEAN PHARMACOPOEIA 10

Pholcodine monohydrate

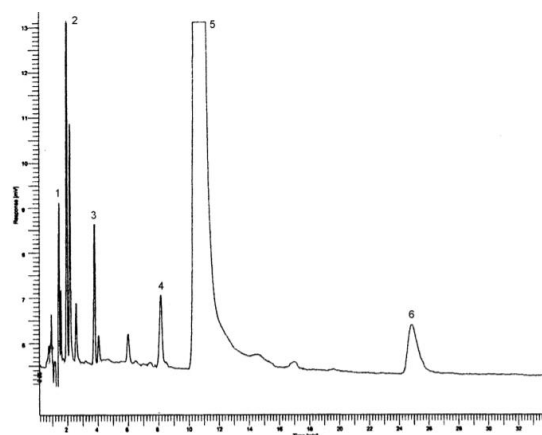


- | | | | | | |
|---------------|---------------|---------------|---------------|---------------|---------------|
| 1. impurity F | 2. impurity C | 3. impurity A | 4. impurity B | 5. pholcodine | 6. impurity D |
|---------------|---------------|---------------|---------------|---------------|---------------|

癸酸睾酮-Testosterone Decanoate

有关物质: Luna C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Testosterone decanoate

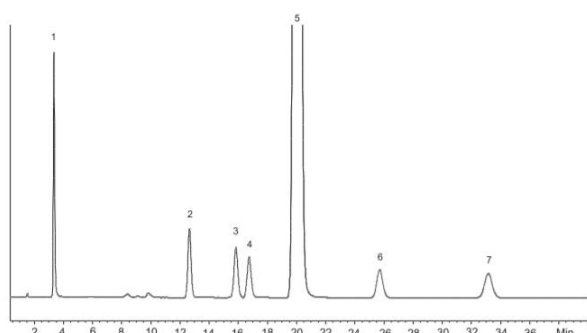


- | | | | |
|---------------|-----------------------|---------------------------|---------------|
| 1. impurity A | 3. impurities C and G | 5. testosterone decanoate | 7. impurity F |
| 2. impurity B | 4. impurity D | 6. impurity E | |

异己酸睾酮-Testosterone Isocaproate

有关物质: Luna C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Testosterone isocaproate

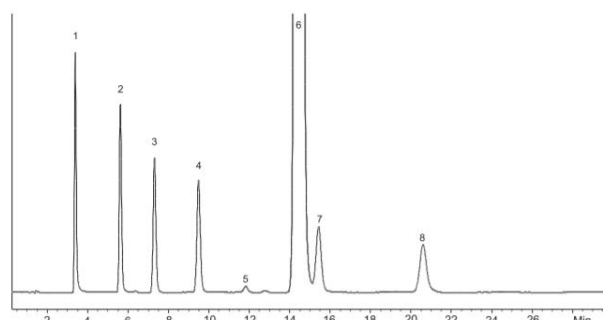


- | | | | |
|---------------|---------------|-----------------------------|---------------|
| 1. impurity A | 3. impurity C | 5. impurity G | 7. impurity E |
| 2. impurity B | 4. impurity D | 6. testosterone isocaproate | 8. impurity F |

庚酸睾酮-Testosterone enantate

有关物质: Synergi MAX-RP (4.6 x150,4μm)

EUROPEAN PHARMACOPOEIA 10

Testosterone enantate

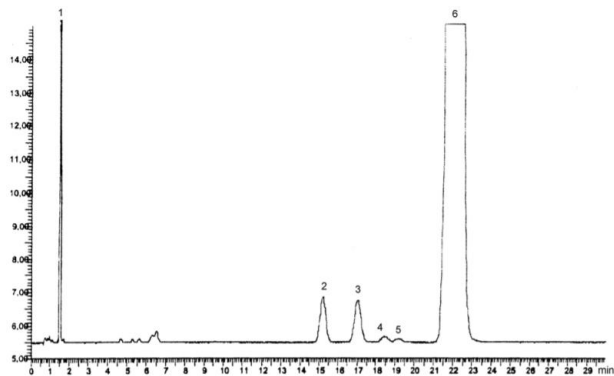


1. impurity D 2. impurity B 3. impurity E 4. impurity F 5. impurity G 6. testosterone enantate

培美曲塞二钠七水合物-Pemetrexed disodium heptahydrate

对映体纯度: IB-Sil C18 (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Pemetrexed disodium heptahydrate

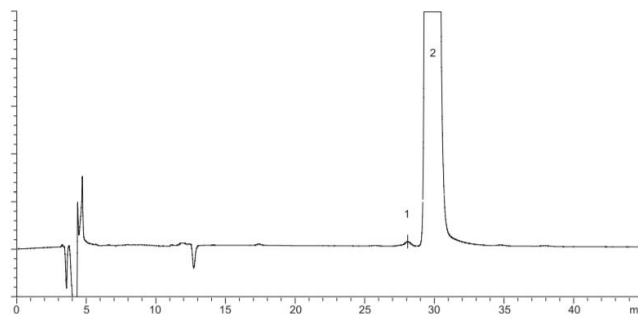


1. impurity E 2. pemetrexed

利塞膦酸钠 2.5 水合物-Risedronate sodium 2.5-hydrate

有关物质: Luna C18(2) (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Risedronate sodium 2.5-hydrate

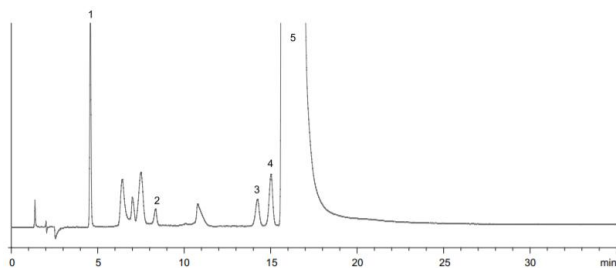


1. impurity D 2. impurity B 3. impurity C 4. impurity E 5. risedronate

硫酸庆大霉素- Gentamicin sulfate

成分和有关物质: Gemini C18 (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Gentamicin sulfate

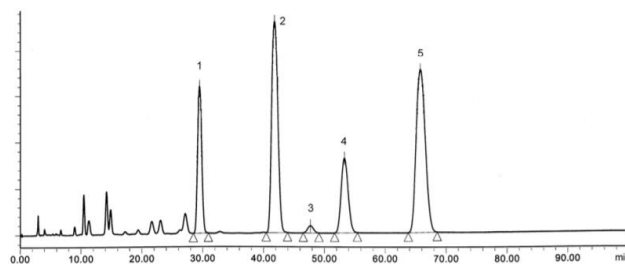


1. gentamicin C1a 3. gentamicin C2b 5. gentamicin C1
2. gentamicin C2 4. gentamicin C2a

硫酸氢氯吡格雷-Clopidogrel

hydrogen sulfate

有关物质: Luna C18 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10 Clopidogrel hydrogen sulfate

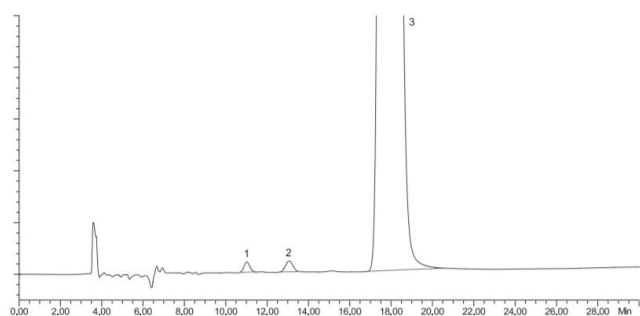


1. impurity C 2. impurity B 3. clopidogrel

吗多明-Molsidomine

有关物质: Luna C18 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Molsidomine

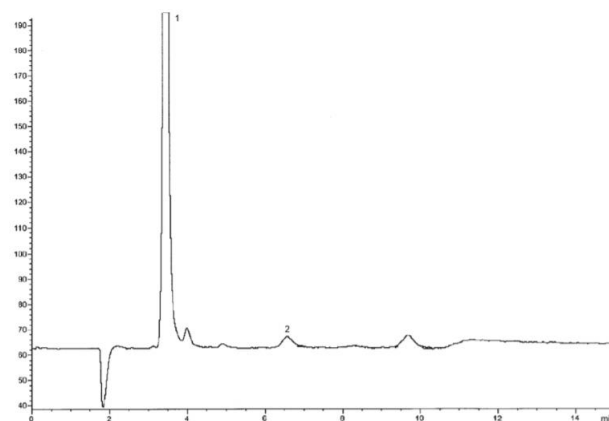


1. molsidomine 2. impurity E

度他雄胺-Dutasteride

有关物质方法 A 和含量: Luna C18
(4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Dutasteride

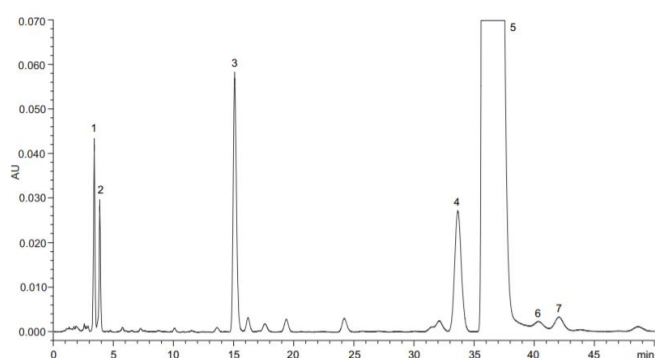


1. impurity A 2. impurity B 3. impurity C 4. impurity E 5. dutasteride 6. impurity F 7. impurity G

米诺地尔-Minoxidil

有关物质: Kinetex XB-C18 (4.6 x100,2.6µm)

EUROPEAN PHARMACOPOEIA 10

Minoxidil

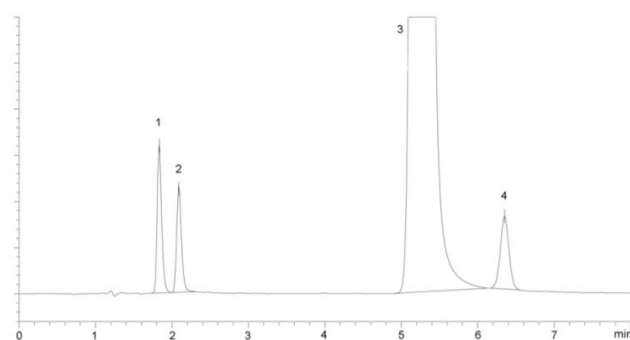


1. impurity A 2. impurity B 3. minoxidil 4. impurity E

2,4-二氯苄醇-2,4-Dichlorobenzyl alcohol

有关物质: Prodigy ODS-3V (4.6 x150,5μm)

EUROPEAN PHARMACOPOEIA 10

2,4-Dichlorobenzyl alcohol

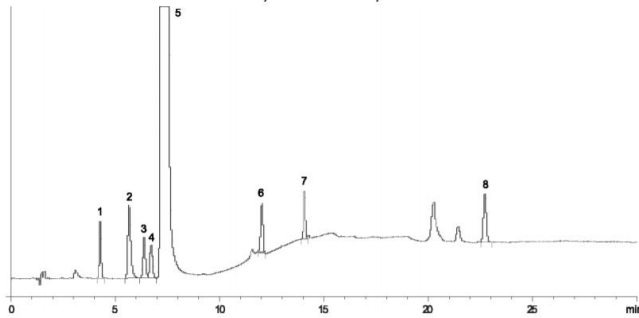


- | | | | |
|---------------|---------------|-------------------------------|---------------|
| 1. impurity B | 3. impurity C | 5. 2,4-dichlorobenzyl alcohol | 7. impurity D |
| 2. impurity E | 4. impurity A | 6. impurity F | 8. impurity G |

卡比多巴-Carbidopa

有关物质: Gemini-NX C18 (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Carbidopa

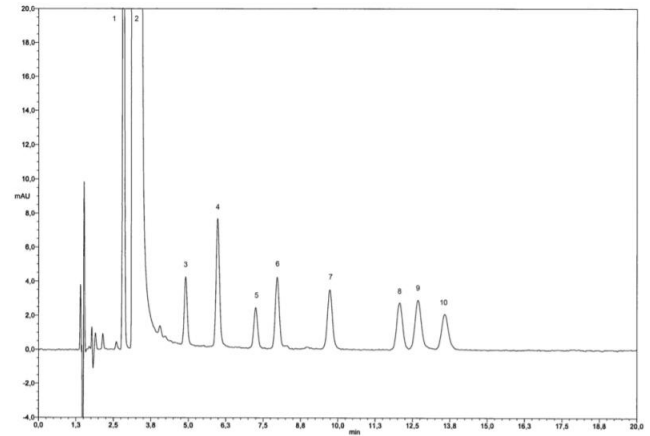


- | | | | | |
|---------------|-------------------|---------------|---------------|----------------|
| 1. impurity A | 3. impurity B | 5. impurity C | 7. impurity G | 9. impurity J |
| 2. carbidopa | 4. impurity D + E | 6. impurity H | 8. impurity I | 10. impurity F |

异麦芽酮糖醇-Isomalt

有关物质: Rezex RCM (7.8 x300)

EUROPEAN PHARMACOPOEIA 10

IsomaltTM

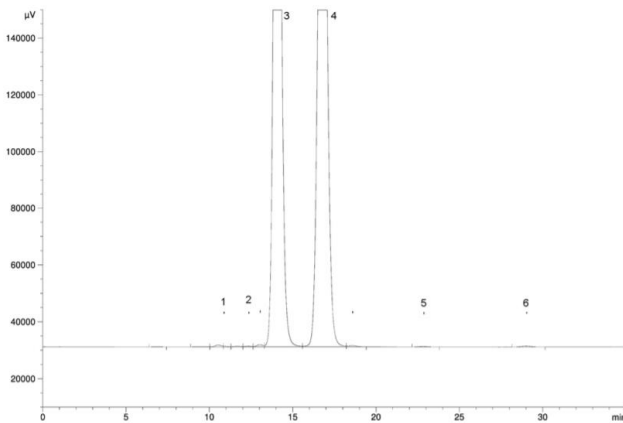

- | | | | | | |
|---------------|---------------|------------|------------|---------------|---------------|
| 1. impurity A | 2. impurity D | 3. 1,1-GPM | 4. 1,6-GPS | 5. impurity B | 6. impurity C |
|---------------|---------------|------------|------------|---------------|---------------|


二水合盐酸纳洛酮-Naloxone hydrochloride dihydrate

hydrochloride dihydrate

杂质 D: Gemini C18 (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Naloxone hydrochloride dihydrate

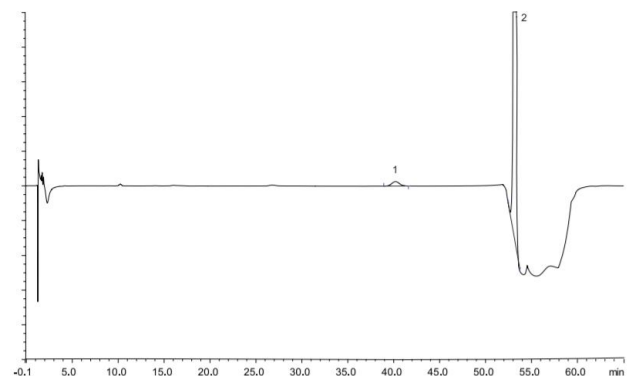


- | | |
|---------------|-------------|
| 1. impurity D | 2. naloxone |
|---------------|-------------|

磺胺二甲氧嘧-Sulfadimethoxine

有关物质: Gemini C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Sulfadimethoxine

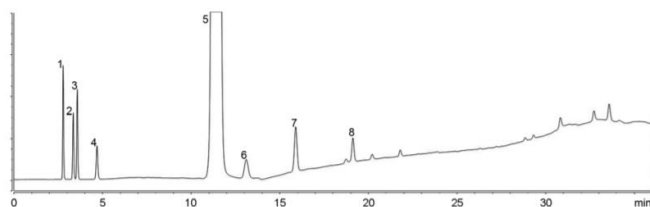

- | | | |
|---------------|---------------------|---------------|
| 1. impurity D | 4. impurity F | 7. impurity B |
| 2. impurity C | 5. sulfadimethoxine | 8. unknown |
| 3. impurity E | 6. impurity A | |


氯己定二葡糖酸盐溶液-

Chlorhexidine digluconate solution

有关物质: Luna C18(2) (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Chlorhexidine digluconate solution

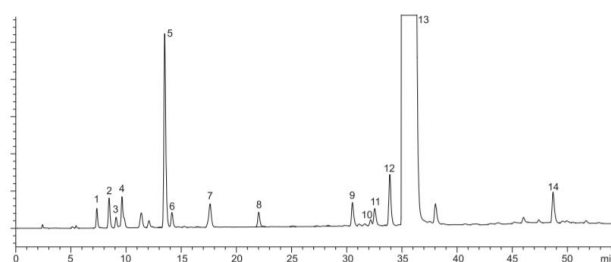


- | | | | | |
|---------------|---------------|---------------|----------------|-------------------|
| 1. impurity E | 4. impurity G | 7. impurity F | 10. impurity O | 13. chlorhexidine |
| 2. impurity L | 5. impurity N | 8. impurity A | 11. impurity I | 14. impurity K |
| 3. impurity Q | 6. impurity B | 9. impurity H | 12. impurity J | |

甲基苯巴比妥-Methylphenobarbital

有关物质: Luna C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Methylphenobarbital

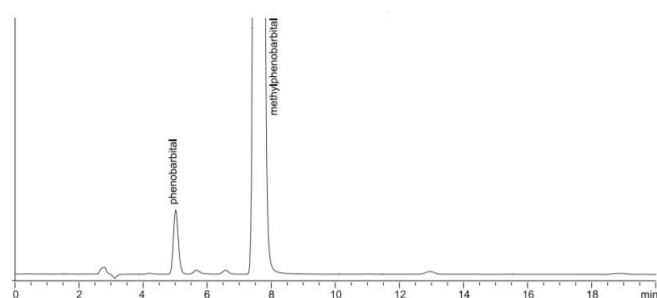


磺胺嘧啶-Sulfadiazine

有关物质: Gemini C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Sulfadiazine

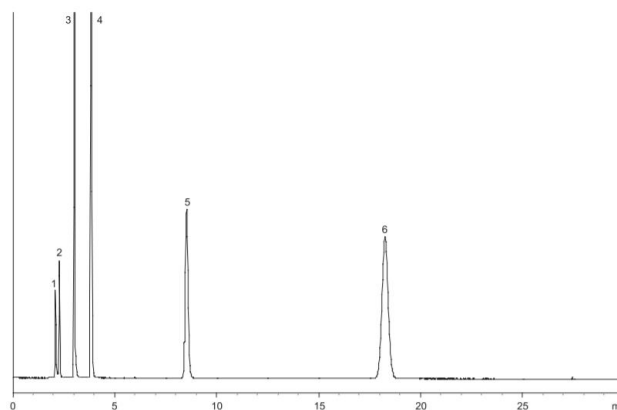


- | | | | | | |
|---------------|---------------|---------------|---------------|-----------------|---------------|
| 1. impurity A | 2. impurity B | 3. impurity C | 4. impurity D | 5. sulfadiazine | 6. impurity E |
|---------------|---------------|---------------|---------------|-----------------|---------------|

甲苯磺丁脲-Tolbutamide

有关物质: Luna C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Tolbutamide

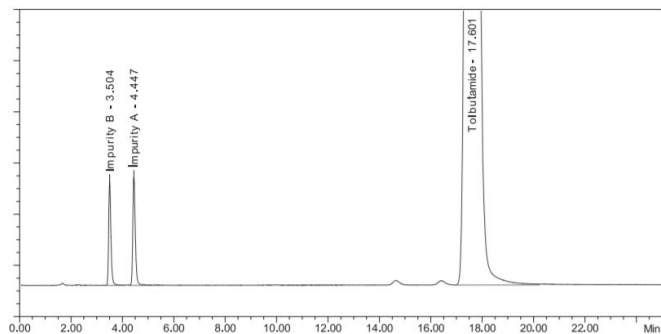


尼泊金甲酯-Methyl

parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Methyl parahydroxybenzoate (1)

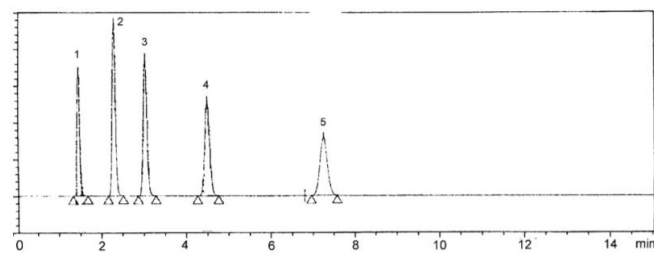

1. impurity A 2. methyl parahydroxybenzoate 3. impurity B 4. impurity C 5. impurity D


尼泊金丙酯-Propyl

parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Propyl parahydroxybenzoate (1)

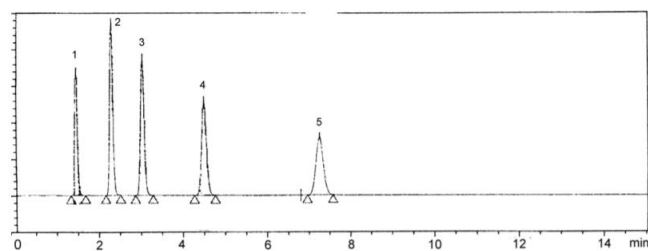


1. impurity A 2. impurity B 3. impurity C 4. propyl parahydroxybenzoate 5. impurity D

烟酸-Nicotinic acid

有关物质: Synergi Hydro-RP (4.6 x250,4µm)

EUROPEAN PHARMACOPOEIA 10

Nicotinic acid

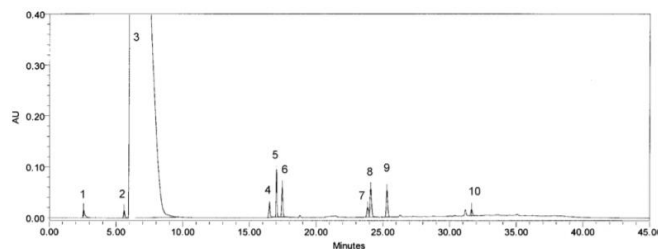

1. impurity D 3. nicotinic acid 5. impurity B 7. impurity G 9. impurity I
2. impurity E 4. impurity A 6. impurity F 8. impurity H 10. impurity C


盐酸布比卡因-Bupivacaine

hydrochloride

杂质 F: Gemini NX-C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Bupivacaine hydrochloride

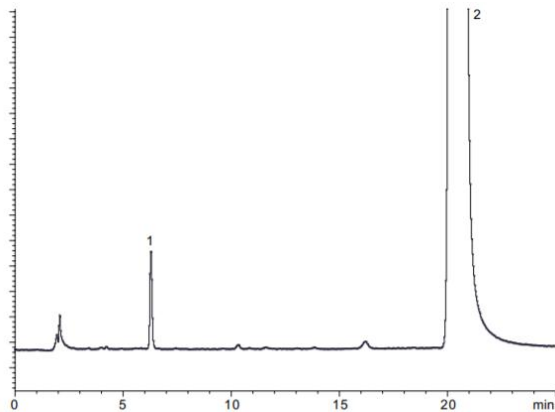

1. impurity F 2. bupivacaine


尼泊金丁酯-Butyl

parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Butyl parahydroxybenzoate (1)

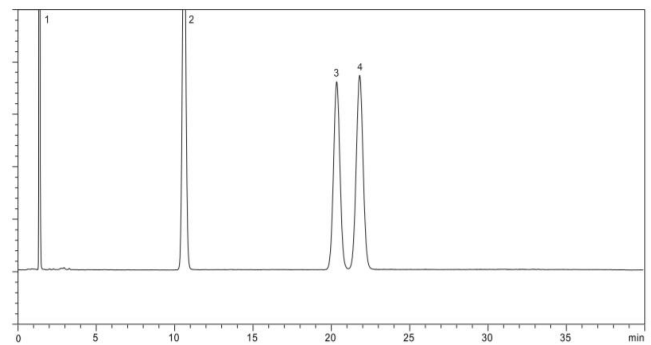

1. impurity A 2. impurity D 3. impurity E 4. butyl parahydroxybenzoate


尼泊金乙酯-Ethyl

parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Ethyl parahydroxybenzoate (1)

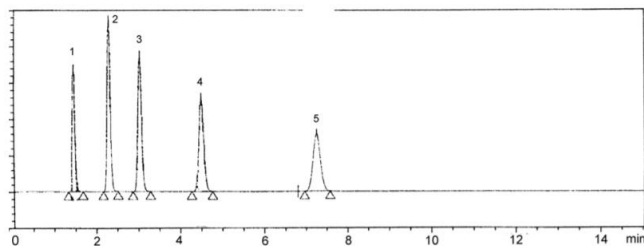

1. impurity A 2. impurity B 3. ethyl parahydroxybenzoate 4. impurity C 5. impurity D


盐酸氟奋乃静-Fluphenazine

dihydrochloride

有关物质: Luna C18 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Fluphenazine dihydrochloride

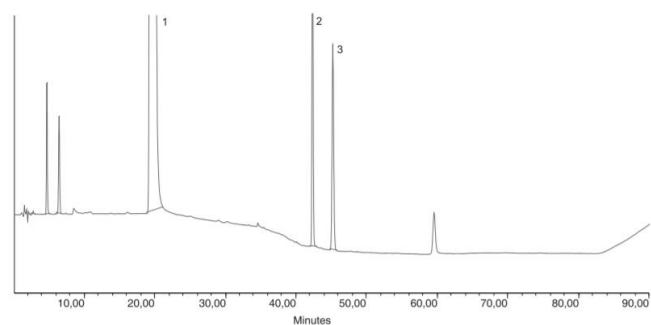


1. fluphenazine 2. impurity D 3. impurity C

尼泊金乙酯钠盐- Sodium ethyl parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Sodium ethyl parahydroxybenzoate

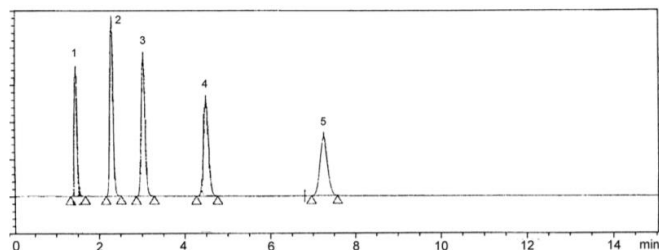


1. impurity A 3. ethyl parahydroxybenzoate 5. impurity D
2. impurity B 4. impurity C

尼泊金甲酯钠盐-Sodium methyl parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 7

Sodium methyl parahydroxybenzoate

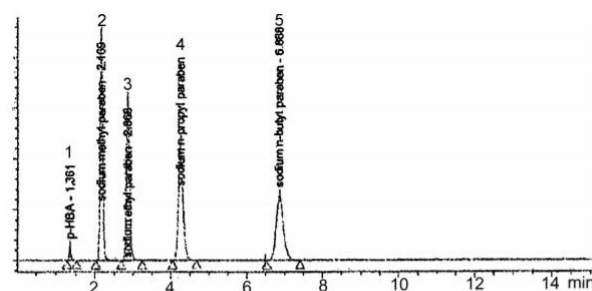


1. impurity A 2. methyl parahydroxybenzoate 3. impurity B 4. impurity C 5. impurity D

尼泊金丙酯钠盐-Sodium propyl parahydroxybenzoate

有关物质: Prodigy ODS-3 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 7

Sodium propyl parahydroxybenzoate

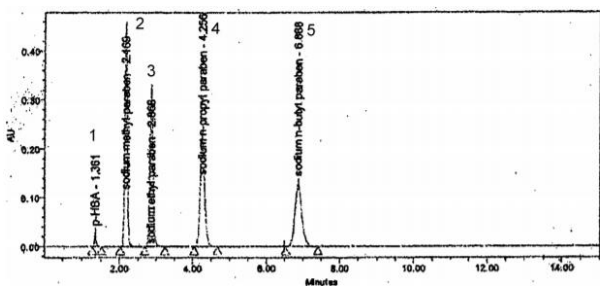


1. impurity A 2. impurity B 3. impurity C 4. propyl parahydroxybenzoate 5. impurity D

醋酸氟卡胺-Flecainide acetate

有关物质: Luna C8 (4.6 x150,5µm)

EUROPEAN PHARMACOPOEIA 10

Flecainide acetate

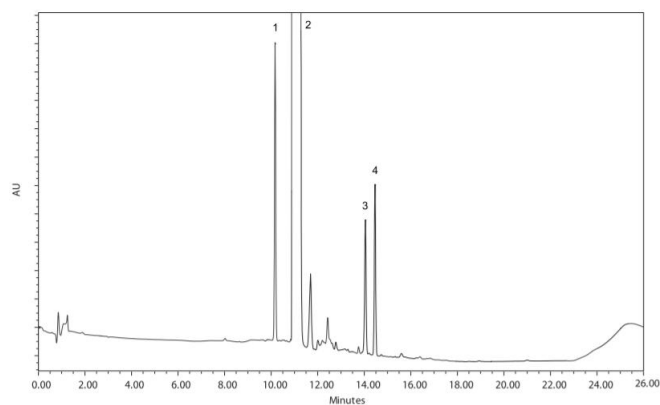


1. impurity C 2. flecainide 3. impurity E 4. impurity D

左卡尼汀- Levocarnitine

有关物质: Sphereclone NH2 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Levocarnitine

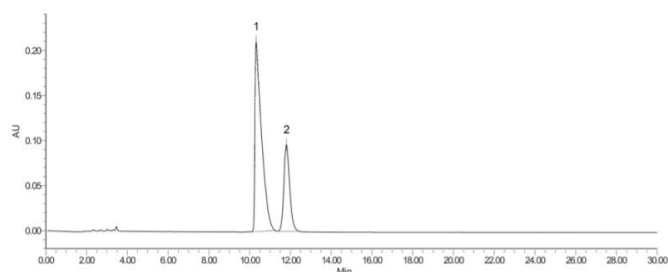

1. levocarnitine


2. impurity A

比卡鲁胺-Bicalutamide

有关物质和含量: Gemini C18 (4.0 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Bicalutamide

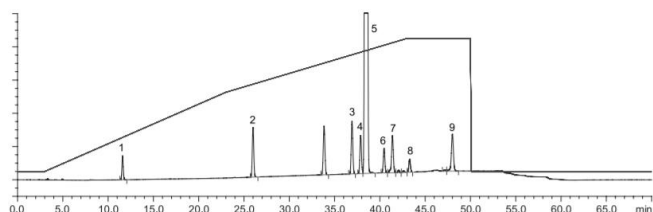

1. impurity M

3. impurity A

5. bicalutamide

7. impurity C

9. impurity J

2. impurity D

4. impurity B

6. impurity K

8. impurity L


氢溴酸西酞普兰-Citalopram

hydrobromide

有关物质: Synergi Hydro-RP (4.6 x250,4µm)

EUROPEAN PHARMACOPOEIA 10

Citalopram hydrobromide

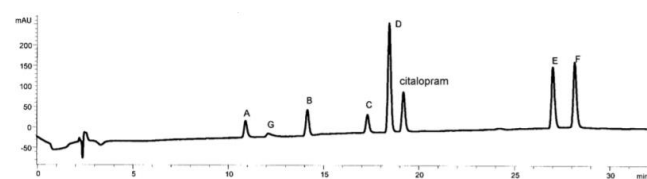


盐酸多佐胺-Dorzolamide

hydrochloride

有关物质: Luna C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Dorzolamide hydrochloride

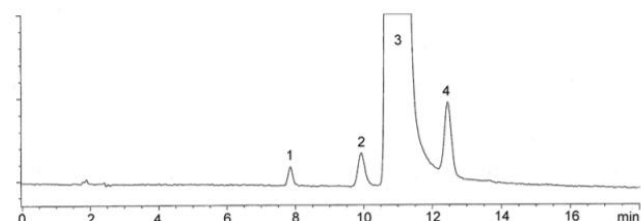

1. impurity D

2. impurity C

3. dorzolamide


4. impurity B

奥卡西平-Oxcarbazepine

有关物质: Luna Phenyl-Hexyl (4.6 x250,5 μ m)

EUROPEAN PHARMACOPOEIA 10

Oxcarbazepine

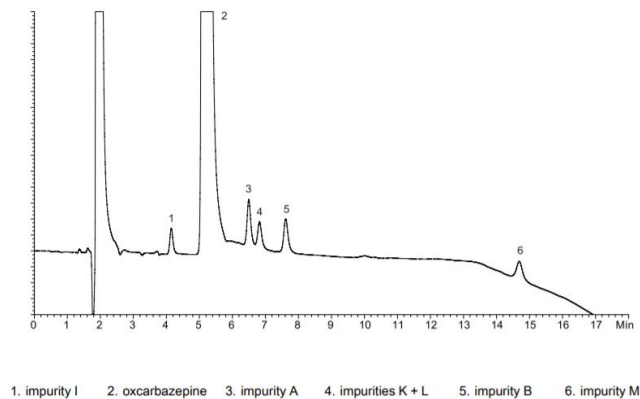


盐酸溴己新-Bromhexine

hydrochloride

有关物质和含量: Kinetex C18 (2.1 x150,2.6 μ m)

EUROPEAN PHARMACOPOEIA 10

Bromhexine hydrochloride

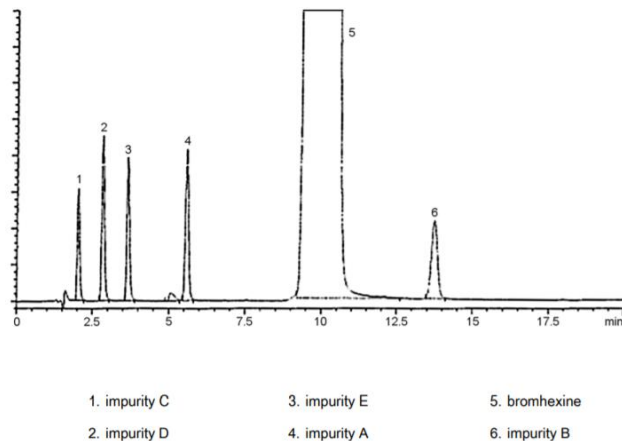


苯巴比妥钠-Phenobarbital sodium

有关物质: Luna C18(2) (4.6 x250,5 μ m)

EUROPEAN PHARMACOPOEIA 10

Phenobarbital sodium

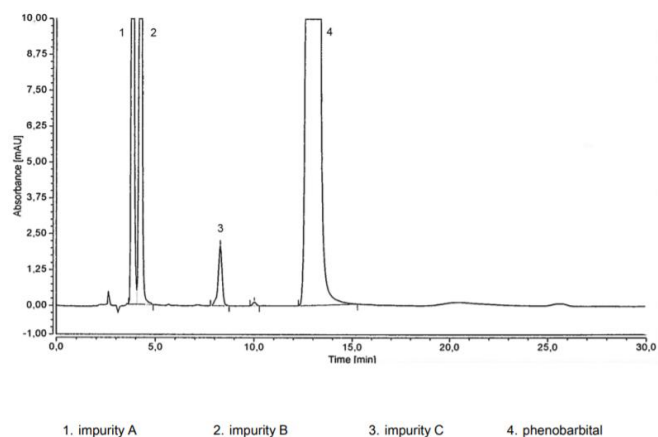


钆布醇一水合物-Gadobutrol

monohydrate

有关物质: Luna Phenyl-Hexyl (4.6 x250,3 μ m)

EUROPEAN PHARMACOPOEIA 10

Gadobutrol monohydrate

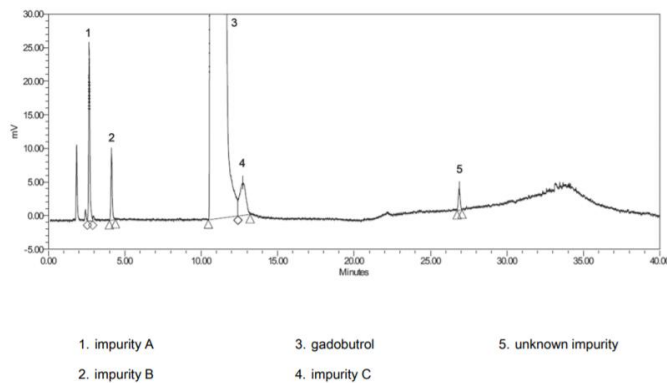


恩替卡韦一水合物-Entecavir

monohydrate

杂质 F: Luna C18 (4.6 x50,5µm)

EUROPEAN PHARMACOPOEIA 10

Entecavir monohydrate

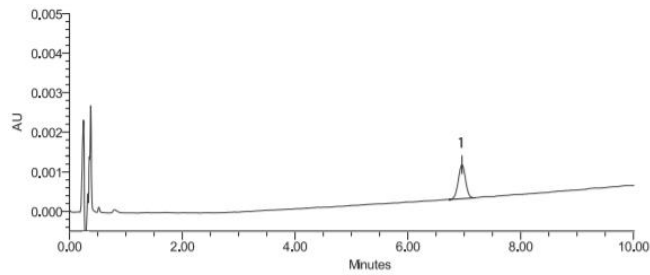

1. impurity F


他卡西醇一水合物- Tacalcitol

monohydrate

有关物质和含量: Luna C18 (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Tacalcitol monohydrate

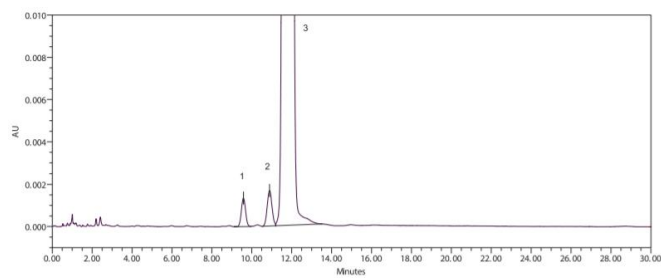

1. pre-tacalcitol

2. impurity A

3. tacalcitol


草酸艾司西酞普兰-Escitalopram

oxalate

有关物质: Luna C18 (2) (4.6 x250,5µm)

EUROPEAN PHARMACOPOEIA 10

Escitalopram oxalate

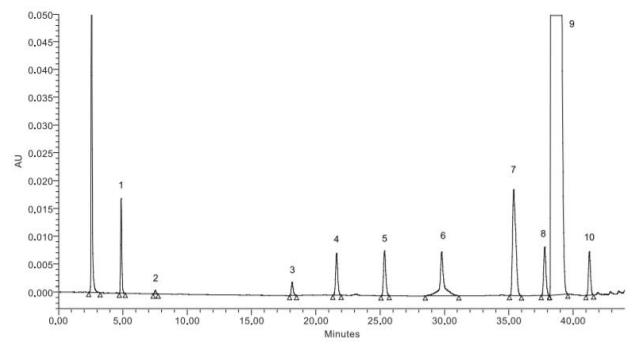

1. impurity F

4. impurity A

7. impurities C+I+L

10. impurity H

2. impurity J

5. impurity E

8. impurity D

3. impurity G

6. impurities B+M


9. escitalopram

甲酸-Formic acid

有关物质: Synergi Hydro RP (4.6 x250,4µm)

EUROPEAN PHARMACOPOEIA 10

Formic acid

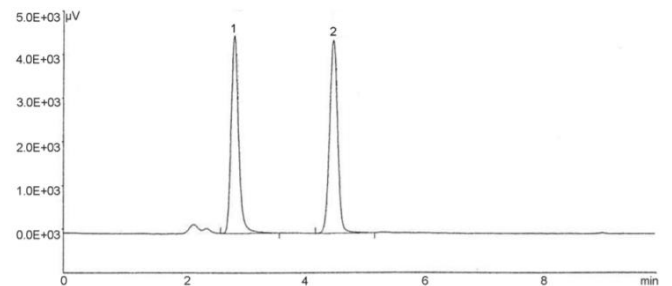

1. formic acid

2. impurity A


他克莫司一水合物-Tacrolimus

monohydrate

有关物质和含量: Luna C18 (2) (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Tacrolimus monohydrate

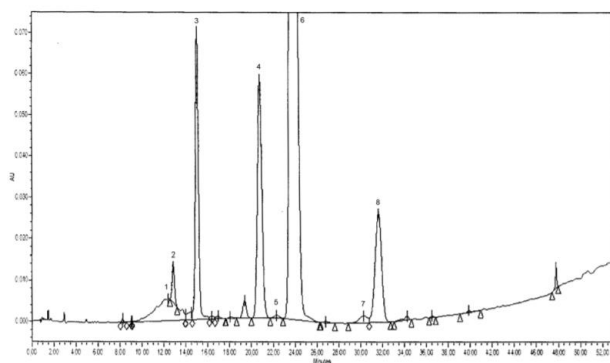


- | | | |
|---------------------------|---------------|---------------|
| 1. tacrolimus compound I | 4. impurity A | 7. impurity D |
| 2. impurity B | 5. impurity C | 8. impurity E |
| 3. tacrolimus compound II | 6. tacrolimus | |

艾司西酞普兰-Escitalopram

有关物质: Luna C18 (2) (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Escitalopram

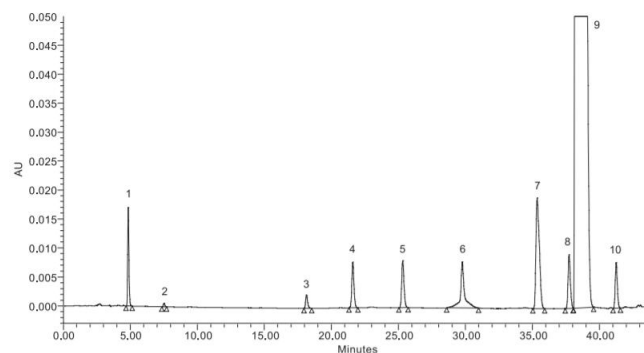


- | | | | |
|---------------|---------------------|-------------------------|----------------|
| 1. impurity F | 4. impurity A | 7. impurities C + I + L | 10. impurity H |
| 2. impurity J | 5. impurity E | 8. impurity D | |
| 3. impurity G | 6. impurities B + M | 9. escitalopram | |

替莫唑胺-Temozolomide

有关物质和含量: Luna C18 (2) (4.6 x150,5μm)

EUROPEAN PHARMACOPOEIA 10

Temozolomide

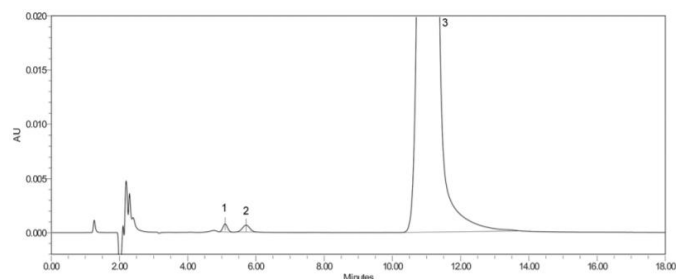


- | | | |
|---------------|---------------|-----------------|
| 1. impurity E | 2. impurity D | 3. temozolomide |
|---------------|---------------|-----------------|

蔗糖-Sucrose, liquid

有关物质: Rezex RCM (7.8 x300)

EUROPEAN PHARMACOPOEIA 10

Sucrose, liquid

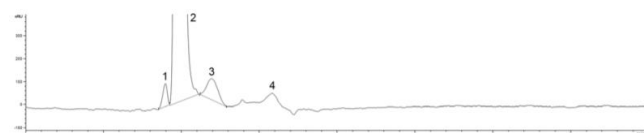

- | | |
|--------------------|---------------|
| 1. impurity A or B | 3. impurity C |
| 2. sucrose | 4. impurity D |


盐酸尼卡地平-Nicardipine

hydrochloride

有关物质: Ultracarb ODS (20) (4.6 x150,5μm)

EUROPEAN PHARMACOPOEIA 10

Nicardipine hydrochloride

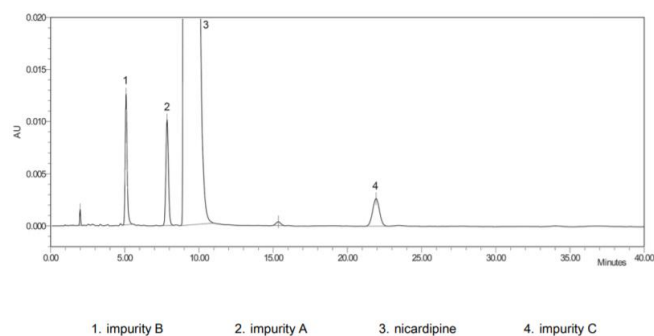


替考拉宁-Teicoplanin

成分和有关物质: Gemini NX-C18
(4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Teicoplanin

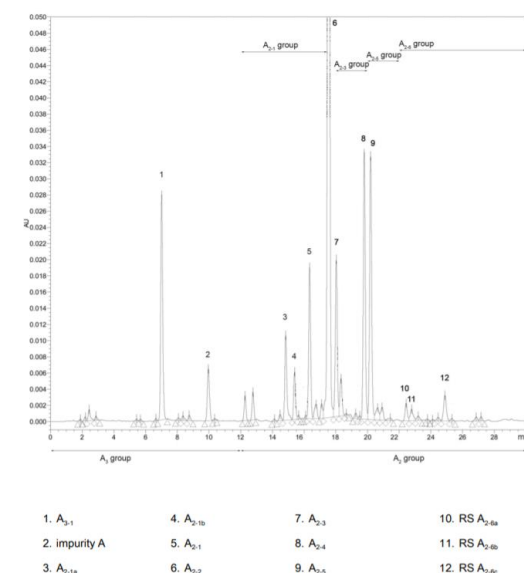


富马酸卢帕他定-Rupatadine

fumarate

有关物质: Gemini NX-C18 (4.6 x150,5μm)

EUROPEAN PHARMACOPOEIA 10

Rupatadine fumarate

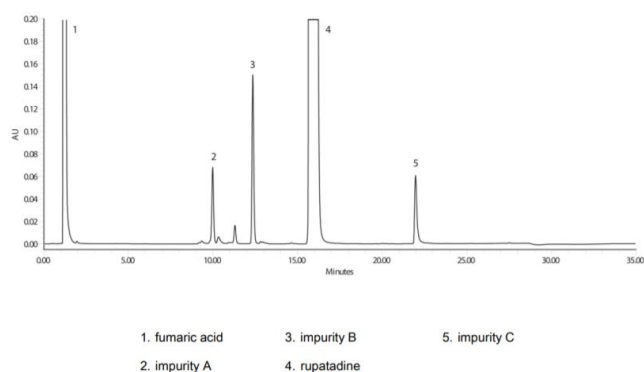


替加环素-Tigecycline

有关物质: Luna C18 (2) (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Tigecycline

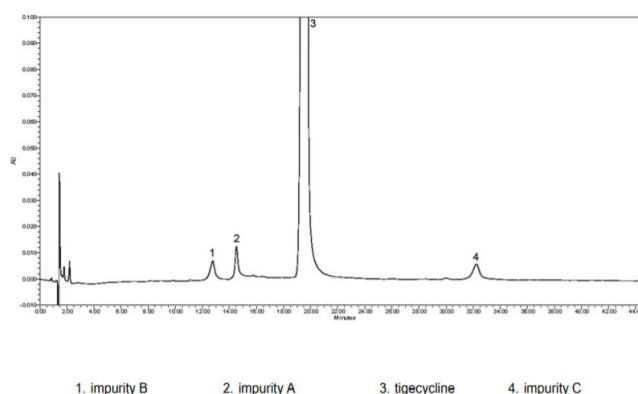


依那西普-Ethosuximide

有关物质: Kinetex EVO C18 (2.1 x100,2.6μm)

EUROPEAN PHARMACOPOEIA 10

Ethosuximide

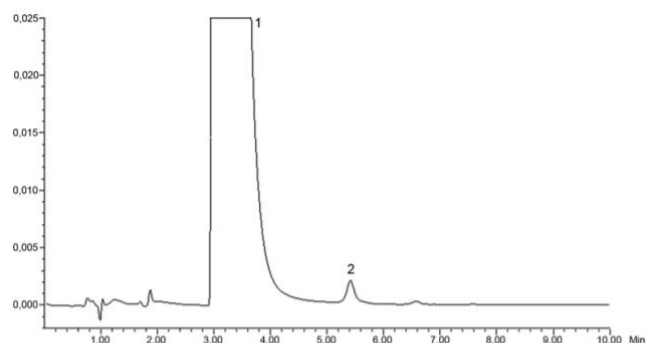

1. ethosuximide


2. impurity A

唑来磷酸一水化合物-Zoledronic acid monohydrate

有关物质: Luna Phenyl-Hexyl (4.6 x250,5μm)

EUROPEAN PHARMACOPOEIA 10

Zoledronic acid monohydrate

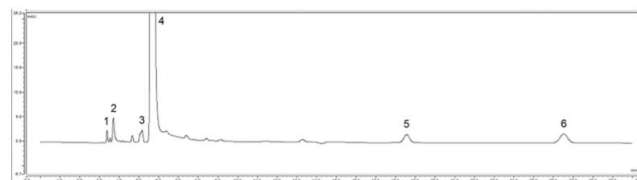

1. nitrate

3. impurity A

5. impurity D

2. impurity B

4. zoledronic acid


6. impurity C

骨化二醇一水合物-Calcifediol monohydrate

有关物质和含量: Luna C8 (4.6 x150,5μm)

EUROPEAN PHARMACOPOEIA 10

Calcifediol monohydrate

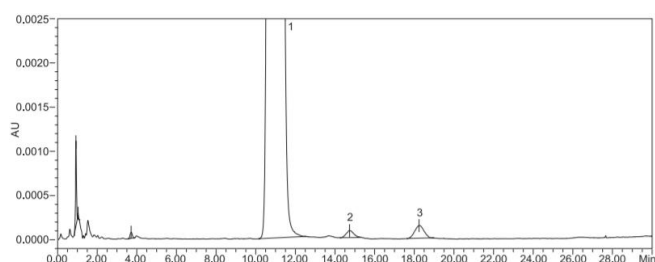

1. calcifediol

2. pre-calcifediol


3. impurity D

盐酸氟哌噻吨-Flupentixol dihydrochloride

有关物质: Gemini-NX C18 (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Flupentixol dihydrochloride

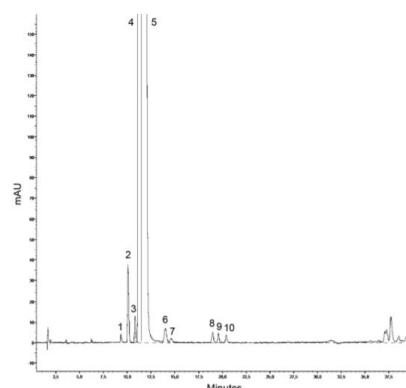

1. impurity A

3. impurities C+I

5. flupentixol (E-isomer)

7. impurity B (isomer 2)

9. impurity E (isomer 1)

2. impurity H

4. flupentixol (Z-isomer)

6. impurity B (isomer 1)

8. impurity G


10. impurity E (isomer 2)

硫酸阿扎那韦-Atazanavir sulfate

有关物质和含量: Luna C18 (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Atazanavir sulfate

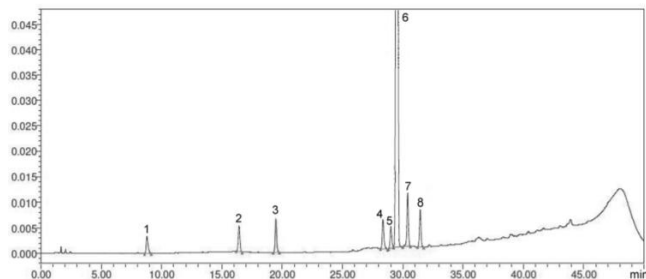

- | | | | |
|---------------|---------------|---------------|---------------|
| 1. impurity A | 3. impurity C | 5. impurity F | 7. impurity G |
| 2. impurity B | 4. impurity E | 6. atazanavir | 8. impurity H |


盐酸多奈哌齐-Donepezil

hydrochloride

有关物质: Luna Phenyl-Hexyl (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Donepezil hydrochloride

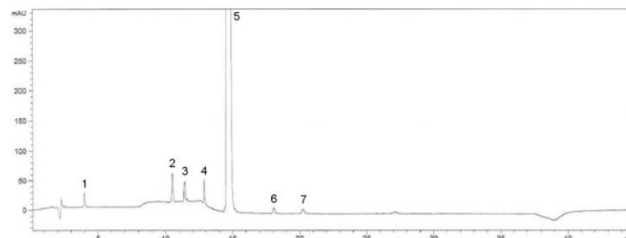

- | | | |
|---------------------|---------------|---------------|
| 1. impurity A | 4. impurity D | 7. impurity G |
| 2. impurity B | 5. donepezil | |
| 3. impurities C + E | 6. impurity F | |


盐酸多奈哌齐一水合物-Donepezil

hydrochloride monohydrate

有关物质: Luna Phenyl-Hexyl (4.6 x150,3μm)

EUROPEAN PHARMACOPOEIA 10

Donepezil hydrochloride monohydrate

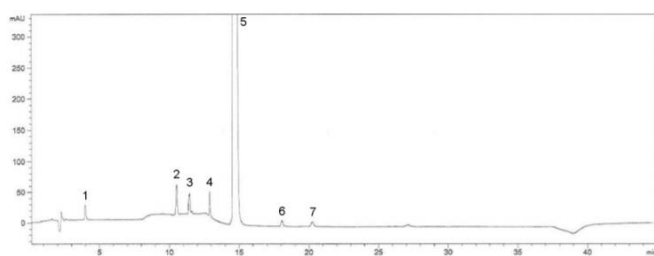

- | | | |
|---------------------|---------------|---------------|
| 1. impurity A | 4. impurity D | 7. impurity G |
| 2. impurity B | 5. donepezil | |
| 3. impurities C + E | 6. impurity F | |


巯嘌呤一水合物-Mercaptopurine

monohydrate

有关物质: Luna C18 (4.6 x100,3μm)

EUROPEAN PHARMACOPOEIA 10

Mercaptopurine monohydrate

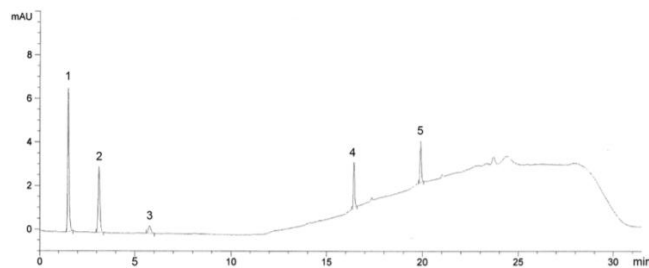


- | | | |
|---------------|-------------------|---------------|
| 1. impurity B | 3. mercaptopurine | 5. impurity D |
| 2. impurity A | 4. impurity C | |

硫酸鱼精蛋白-Protamine sulfate

有关物质：Aeris Widepore XB-C18
(4.6 x150,3.6μm)

EUROPEAN PHARMACOPOEIA 10

Protamine sulfate

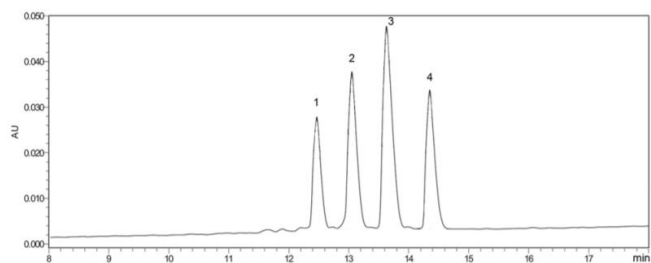

1. protamine A

2. protamine B

3. protamine C

4. protamine D