

# APPLICATIONS

# Fast Screening Method for 86 Forensic Analytes in Human Urine by LC-MS/MS

SCIEX, 1201 Radio Road, Redwood City, CA 94065 USA Phenomenex, Inc., 411 Madrid Ave., Torrance, CA 90501 USA

## **Overview**

This application demonstrates a rapid screening method of 86 compounds in human urine using a Luna® C18(2) LC column and a SCIEX® QTRAP® 4500 MS/MS system. The combination of liquid chromatography coupled with tandem mass spectrometry (LC-MS/MS) allows for simultaneous screening of across multiple compound classes

## **Sample Preparation**

- 100  $\mu$ L urine sample was mixed with 25  $\mu$ L IMCS RapidHydrolysis Buffer, 20  $\mu$ L IMCSzyme and 10  $\mu$ L IS. Hydrolysis time was typically between 30 and 60 min at 55°C.
- After hydrolysis was complete, 0.8 mL of diluent was added to the mixture.
- The mixture was then centrifuged at 21,000 g for 15 min.
- The supernatant was transferred to a glass vial with insert for analysis by LC-MS/MS.

## **LC-MS/MS Conditions**

Column: Luna® 5 μm C18(2)

**Dimension:** 50 x 2.0 mm **Part No.:** 00B-4252-B0

Mobile Phase: A: 5 mM Ammonium formate in water

B: 0.1 % Formic acid in methanol

Gradient: Time (min) %B

0.0 5 1.8 95

Flow Rate: 0.2 - 1.0 mL/min

**Temperature:** Ambient **Injection:** 5 μL

**Detection:** MS/MS (SCIEX® QTRAP® 4500) **Sample:** See Table 1 for list of 86

compounds

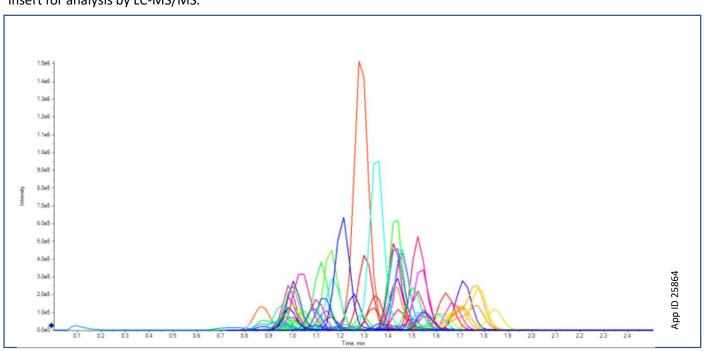




Table 1. 86 Compounds and Internal Standards

Compounds	Compounds	Compounds	Internal Standards
6-MAM	Gabapentin	Naltrexone	Amphetamine-d5
7-Aminoclonazepam	Hydrocodone	N-desmethyltapentadol	Fentanyl-d5
7-Hydroxymitragynine	Hydromorphone	Norbuprenorphine	JWH 018 4-OH pentyl-d5
Acetyl Fentanyl	Imipramine	Norcodeine	Methadone-d3
Alpha-Hydroxyalprazolam	JWH-018 4-OH pentyl	Nordiazepam	Methamphetamine-d5
Alpha-Hydroxymidazolam	JWH-018 pentanoic acid	Norfentanyl	Morphine-d6
Alpha-Hydroxytriazolam	JWH-019 6-OH hexyl	Norhydrocodone	
Alpha-PPP	JWH-073 3-OH butyl	Normeperidine	
Alpha-PVP	JWH-073 butanoic acid	Noroxycodone	_
Alprazolam	JWH-081 5-OH pentyl	Norpropoxyphene	
AM-2201 4-OH pentyl	JWH-122 5-OH pentyl	Nortriptyline	_
Amitriptyline	JWH-210 5-OH pentyl	O-Desmethyltramadol	
Amphetamine	JWH-250 4-OH pentyl	Oxazepam	
Benzoylecgonine	Lorazepam	Oxycodone	_
Buphedrone	MDA	Oxymorphone	
Buprenorphine	MDEA	PCP	
Carisoprodol	MDMA	Pregabalin	-
Clomipramine	MDPV	Propoxyphene	-
Codeine	Meperidine	Protriptyline	
Cotinine	Mephedrone	RCS4-4-OH-pentyl	
Cyclobenzaprine	Meprobamate	Ritalinic Acid	
Desalkylflurazepam	Methadone	Sufentanil	_
Desipramine	Methamphetamine	Tapentadol	
Desmethyldoxepin	Methedrone	Temazepam	
Dextromethorphan	Methylone	Tramadol	
Diazepam	Methylphenidate	Zolpidem	
Dihydrocodeine	Midazolam		
Doxepin	Mitragynine		
EDDP	Morphine		
Fentanyl	Naioxone		

## Key Features of Luna® LC Media Paired with QTRAP® Technology

- Rapid, robust approach allows direct detection of 86 forensic analytes in human urine with simple sample preparation and fast chromatography.
- Scheduled MRM™ Algorithm enables the use of fast chromatography without loss of data integrity, maintaining sufficient data points across the peak for reliable quantification and higher throughput
- Targeted MRM workflow allows sensitive detection with high selectivity, reduced matrix effects and higher confidence in results.
- Additional QTRAP technology allows for enhanced scan functions that further improve confidence in results, by generating high sensitivity MS/MS spectra to enhance confidence of identification of analytes, enabling spectral identification and library searching.

## phenomenex ...breaking with tradition<sup>™</sup>

## PPLICATIONS

## Need a different column size or sample preparation format?

No problem! We have a majority of our available dimensions up on www.phenomenex.com, but if you can't find what you need right away, our super helpful Technical Specialists can guide you to the solution via our online chat portal www.phenomenex.com/ChatNow.

Australia t: +61 (0)2-9428-6444 auinfo@phenomenex.com

### Austria

t: +43 (0)1-319-1301 anfrage@phenomenex.com

**Belgium** t: +32 (0)2 503 4015 (French) t: +32 (0)2 511 8666 (Dutch) beinfo@phenomenex.com

#### Canada

t: +1 (800) 543-3681 info@phenomenex.com

China t: +86 400-606-8099 cninfo@phenomenex.com

### Denmark

t: +45 4824 8048 nordicinfo@phenomenex.com

## **Finland**

t: +358 (0)9 4789 0063 nordicinfo@phenomenex.com

## France

t: +33 (0)1 30 09 21 10 franceinfo@phenomenex.com

Germany t: +49 (0)6021-58830-0 anfrage@phenomenex.com

## India

t: +91 (0)40-3012 2400 indiainfo@phenomenex.com

#### Ireland

t: +353 (0)1 247 5405 eireinfo@phenomenex.com

Italy t: +39 051 6327511 italiainfo@phenomenex.com

## Luxembourg

t: +31 (0)30-2418700 nlinfo@phenomenex.com

#### Mexico

t: 01-800-844-5226 tecnicomx@phenomenex.com

#### The Netherlands

t: +31 (0)30-2418700 nlinfo@phenomenex.com

New Zealand t: +64 (0)9-4780951 nzinfo@phenomenex.com

Norway t: +47 810 02 005 nordicinfo@phenomenex.com

t: +48 (12) 881 0121 pl-info@phenomenex.com

Portugal t: +351 221 450 488 ptinfo@phenomenex.com

**Singapore** t: +65 800-852-3944 sginfo@phenomenex.com

**Spain** t: +34 91-413-8613 espinfo@phenomenex.com

#### Sweden

t: +46 (0)8 611 6950 nordicinfo@phenomenex.com

## Switzerland

t: +41 (0)61 692 20 20 swissinfo@phenomenex.com

#### Taiwan

t: +886 (0) 0801-49-1246 twinfo@phenomenex.com

United Kingdom t: +44 (0)1625-501367 ukinfo@phenomenex.com

USA t: +1 (310) 212-0555 info@phenomenex.com

## All other countries/regions

Corporate Office USA t: +1 (310) 212-0555 info@phenomenex.com

## www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country/region, contact Phenomenex USA, International Department at international@phenomenex.com

Subject to Phenomenex Standard Terms and Conditions, which may be viewed at www.phenomenex.com/TermsAndConditions.

## **Trademarks**

Luna is a registered trademark Phenomenex. SCIEX and QTRAP are registered trademarks and Scheduled MRM is a trademark of AB SCIEX Pte. Ltd. AB SCIEX m is being used under license.

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures. © 2020 Phenomenex, Inc. All rights reserved.